

## KB00022 - ERROR - Sample Master® v8.5 Sample Login Displays ODBC Error 260 When Processing Prelog2Login

### SYMPTOMS:

The Sample Master® V85 Sample Login function displays a message like the one below when a user attempts to process a Prelog to a Login.

ODBC Error 260

Disallowed implicit conversion from data type datetime to data type bit, table 'SMV8.SMSU.vw\_Orders\_Prelog2Login', column 'ReceiveFlag'. Use CONVERT function to run this query.

An error occurred running

SMV8.SMSU.sp\_Orders\_Prelog2Login '2010-07-01 09:39:17

### CAUSE:

This error is caused by a bug in the SMSU.sp\_Orders\_Prelog2Login stored procedure in SQL Server.

### CORRECTIVE ACTION:

The code error is corrected in SMV85\_BE\_SQL\_SP02.EXE BE component Service Pack. Application of all three components of Service Pack SP02 will eliminate the error.

## Applies To:

- Sample Master® v8.5 Pro
- Sample Master® v8.5

### FOR MORE INFO

Contact ATL Support:

- Call: 1-800-565-LIMS (5467)
- Email: [support@atlab.com](mailto:support@atlab.com)
- Monday thru Friday 8 am – 5 pm ET

## Accelerated Technology Laboratories Inc.

Founded in 1994, Accelerated Technology Laboratories, Inc., launched the first commercially available Windows-based Laboratory Information Management System (LIMS). More than two decades and several hundred installations later, Sample Master® is a market-leading solution with thousands of users around the world. ATL is recognized for our strong commitment to quality, investment in research and development, and team of professionals. We hire experts with impeccable academic credentials in the fields of chemistry, engineering, water and wastewater, microbiology, medical technology, toxicology, biotechnology, laboratory management, validation, software development, computer science, engineering, and business. Several hold MS and Ph.D. degrees. ATL professionals are highly sought-after speakers and thought leaders. We are actively involved in the scientific community. Our affiliations and partnerships allow us to deliver the most current leading-edge technology to our users. ATL also offers TITAN LIMS, enhancement products, support, and consulting services worldwide.

For more information: [www.atlab.com](http://www.atlab.com)



496 Holly Grove School Rd  
West End, NC 27376

toll-free: 800.565.LIMS (5467)  
phone: 910.673.8165  
fax: 910.673.8166

[www.atlab.com](http://www.atlab.com)  
[info@atlab.com](mailto:info@atlab.com)