


KB00030 - ERROR - ODBC Error 257 Implicit Conversion of data type DateTime to Bit is not allowed

SYMPTOMS:

In Sample Master® V85 a message like the one below is displayed when a user attempts to transfer a Prelog order to a logged order.

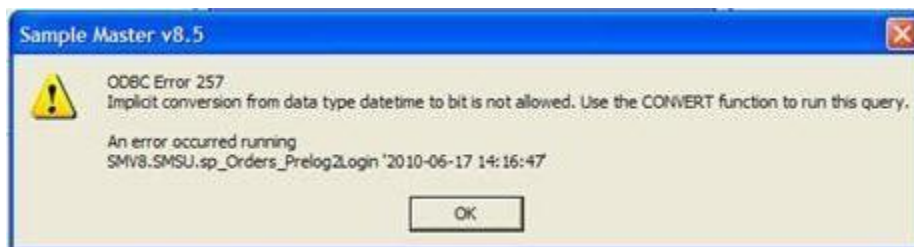

CAUSE:

This error is caused by a bug in the application

CORRECTIVE ACTION:

The error has been resolved in service pack 02. Application of the three components of this service pack will eliminate this error.

Applies To:

- Sample Master® v8.5

FOR MORE INFO

Contact ATL Support:

- Call: 1-800-565-LIMS (5467)
- Email: support@atlab.com
- Monday thru Friday 8 am – 5 pm ET

Accelerated Technology Laboratories Inc.

Founded in 1994, Accelerated Technology Laboratories, Inc., launched the first commercially available Windows-based Laboratory Information Management System (LIMS). More than two decades and several hundred installations later, Sample Master® is a market-leading solution with thousands of users around the world. ATL is recognized for our strong commitment to quality, investment in research and development, and team of professionals. We hire experts with impeccable academic credentials in the fields of chemistry, engineering, water and wastewater, microbiology, medical technology, toxicology, biotechnology, laboratory management, validation, software development, computer science, engineering, and business. Several hold MS and Ph.D. degrees. ATL professionals are highly sought-after speakers and thought leaders. We are actively involved in the scientific community. Our affiliations and partnerships allow us to deliver the most current leading-edge technology to our users. ATL also offers TITAN LIMS, enhancement products, support, and consulting services worldwide.

For more information: www.atlab.com



496 Holly Grove School Rd
West End, NC 27376

toll-free: 800.565.LIMS (5467)
phone: 910.673.8165
fax: 910.673.8166

www.atlab.com
info@atlab.com